

Particle Line Connection

If you have not read the **General Plumbing Instructions** for replacing plumbing parts on the KinExA® please do so now before continuing.

The following **Parts and/or Tools** are required to complete this installation:

[1] Particle Line Connection (344270)

[1] 5/16" Crowfoot Wrench (043050)

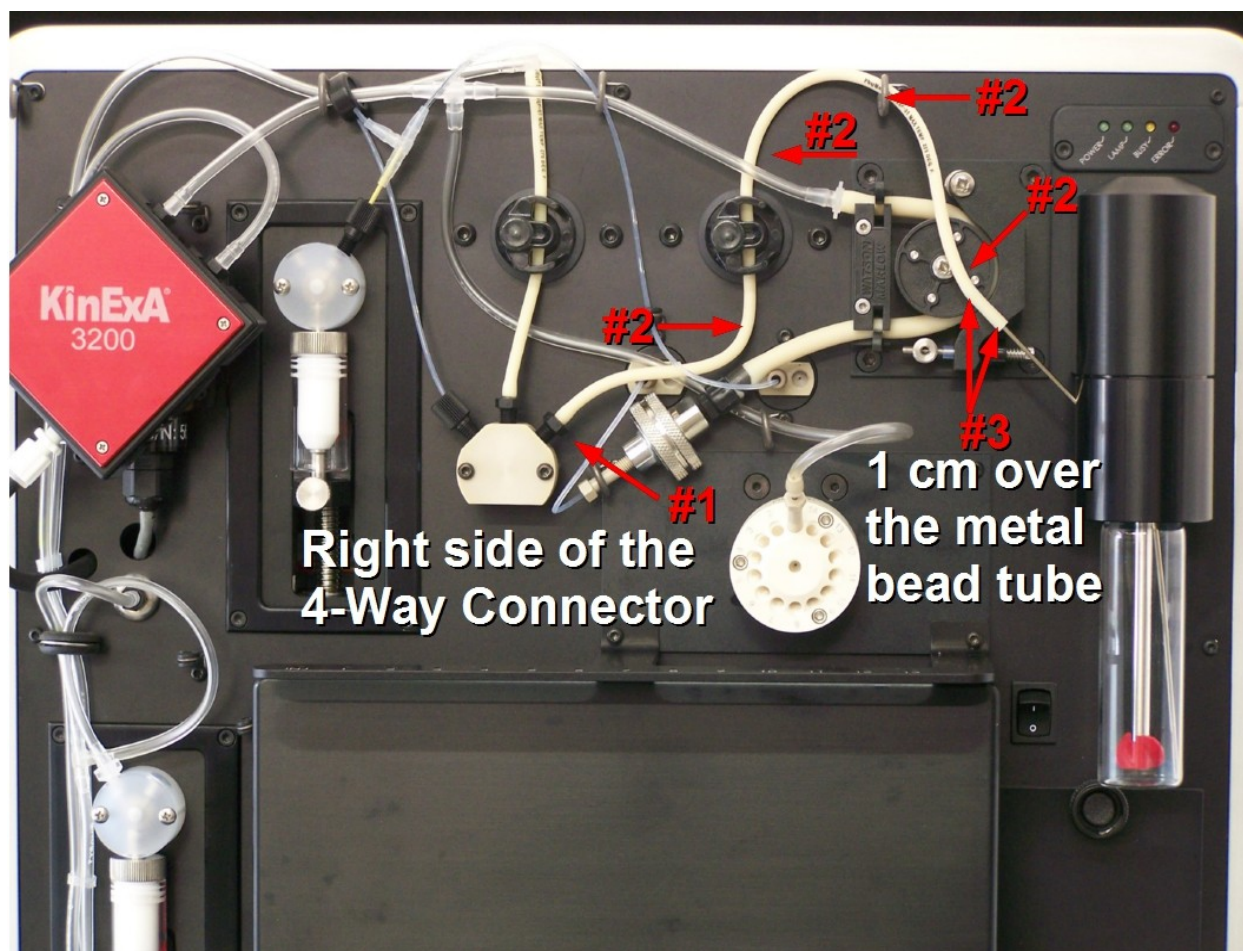


Figure 1: Particle Line Connection routing diagram.

- Screw the *Hose Barb Connector* into the right side of the *4-Way Connector* using the *5/16" Crowfoot Wrench* (043050) (**Figure 1, #1**).
Note: Always hand tighten the *Hose Barb Adapters* followed by a 1/4 turn to insure proper tightness.
- Route the other end over the *Backflush Tubing* above the left *Solenoid Valve*, under the *Injection T Tubing*, and into the right *Pinch Valve*. Continue through the top-right *Tubing Support* and 2 cm over the metal bead tube (**Figure 1, #2**).
- Gently and evenly pull the tubing back to **1 cm** to assure proper alignment (**Figure 1, #3**).